SONO SWISS[®]

High Quality Ultrasonic Systems

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The Spirit of the brand is swissmade

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The new ultrasonic units for finest cleaning applications with aqueous cleaning solutions in the laboratory, in the medical sector, in watchmakers' and jewellers' workshops and in the industrial sector.

We want you to be able to "clean without any residues". The modern ultrasonic cleaning technology of Sonoswiss meets the high demands of today and exceeds all conventional ultrasonic cleaning methods. It is suitable for the cleaning process itself and for further rinsing baths. Sonoswiss Ultrasound with powerful piezo-transducer systems and microprocessor-controlled generators.

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Easy handling of time, temperature and special functions. Clear indication of set and actual values plus visualization by diodes of the special functions.

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*Useful and practical accessories.
The cover can also serve as drip-off collecting basin for the wet basket.*

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*Connection for drain duct at the back, therefore the workplace can be chosen with or without a wash basin.
Comfortable operating elements on the unit side.*

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The housing, made completely of stainless steel, allows hygienic cleaning and disinfection (if required) of the unit surfaces.

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*Protecting rim to keep the operating panel free of cleaning medium.
The operating elements themselves are protected against splash water – in addition any spilled water is guided past the operating panel.*

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For the cleaning of: glassware and complicated glass/ceramics instruments and parts, metal and plastic.

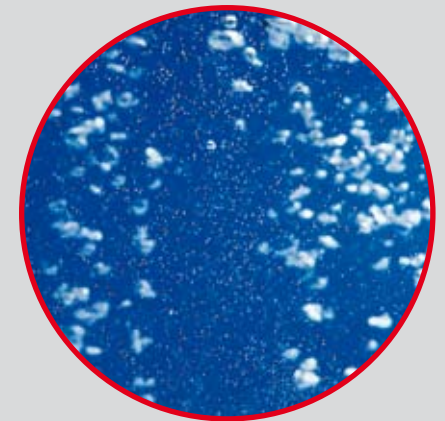
Removal of: chemical residues, coking and soot, baked residues of chemical reactions.

Frequent laboratory applications:

Sonoswiss ultrasonic cleaning units are indispensable helpers for everyday laboratory procedures. Glassware and instruments treated with a suitable cleaning medium will be as clean as new with no traces of contamination left behind, so results of analyses cannot be falsified.

In addition to cleaning jobs, the units are used for the degassing of samples, for decarbonating and for the degassing of HPLC solutions. Furthermore, the units are perfect for removing hardly soluble substances, and for dispersing and homogenizing.

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Degas: samples and HPLC solutions must be free of gas bubbles as far as possible. Quick degassing without the formation of bubbles.

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For the cleaning of: *used surgical instruments in the hospital, in the medical and dentist's practice.*

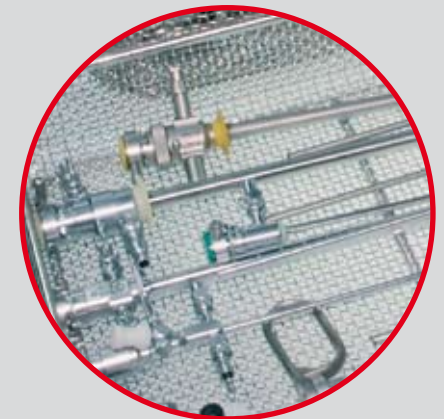
Removal of: *organic contaminants, sterilizing residues, silicate stains from rinsing media and other substances.*

In the dental laboratory impression spoons, prostheses and dental drills are reliably cleaned without any residues.

Sonoswiss units are a fix constituent in the hygienic chains of hospitals and medical practices.

Due to our high demand on the cleaning results our units guarantee the full use of the material service life of the instruments and protect the full range of their functions.

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Sweep: *Instruments must be cleaned reliably no matter where they are placed in the bath. The Sweep function guarantees an even sound field distribution.*

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For the cleaning of: watches, worn or repaired jewellery, stored pieces and jewellery in the jeweller's workshop.

Removal of: oxide layers (layers forming during prolonged periods of storing or after soldering), polishing pastes (after repair or in the production process) and grease and traces of make-up on worn jewellery. Our units are particularly suitable for the cleaning of pieces made of gold, platinum/iridium and silver, also for hard precious stones such as diamonds and sapphires. Caution with soft stones and corals! These could be destroyed or delustered. Watchmakers use our units for the cleaning of parts of clocks and pendulums. Resinified oils and oxide layers are removed. For small watches we recommend to use our units only to clean the metal straps.

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Heating: grease and wax are removed in aqueous cleaning media only when heated above their fusion temperature. Cleaning solutions must be heated.

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For the cleaning of: *productional parts in industrial and handicraft shops. Special pieces that require further processing after the cleaning, e.g. with galvanic plating, titanium nitride (TiN) coating must first be removed without residue and must be conditioned correctly. Parts such as micro-mechanics, pieces with bores or otherwise barely accessible grooves or apertures must be deaerated after immersion.*

Removal of e.g.: *lapping and polishing media, grease, oil, wax, scale. The Sonoswiss units are well suitable for the tough application in the industrial manufacturing processes. The splash water protection rim protects the operating elements from liquid and contaminations. The cover (optional accessory) is made of polypropylene and very resistant against aqueous cleaning media.*

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Boost: *the ultrasound regime is modified which increases the cleaning effect for tenacious contaminations.*

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*We offer a large range of useful and practical accessories for our ultrasonic cleaning units. Please ask for the list of accessories!
 You will obtain perfect cleaning results by the use of special aqueous cleaning concentrates, either liquid or pulverized. Please ask for our list of recommended cleaning media!*

SONO SWISS

Sonoswiss AG
 Sonnenstr. 417
 CH-8262 Ramsen
 Telefon +41 (52) 742 80 10
 Telefax +41 (52) 742 80 18
 info@sonoswiss.ch
 www.sonoswiss.ch